

There are two DataPlate® digital hotplates [Fig 1] in the Microfabrication cleanroom located next to the SCS Spincoater.

Typical Operating Procedure

1. Switch on main power. [Fig 2]
2. After a few seconds, current plate temperature will display.

Set-Point

1. Press SET button. [Fig 3]
2. Press PLATE TEMP button.
3. Input set-point temperature.
 - Note:** Due to age the 1 and 2 button faces are worn.
 - PLATE TEMP = 1 and PROBE TEMP = 2.
4. Press SET button to accept inputted set-point temperature.
 - Note:** The set-point temperature and plate temperature will display sequentially on digital display.
5. To change plate temperature, repeat steps 1-4.

CAUTION: Always turn off hotplate by switching off main power [Fig 2] when not in use. Never leave hotplate unattended when in use.



Fig 1

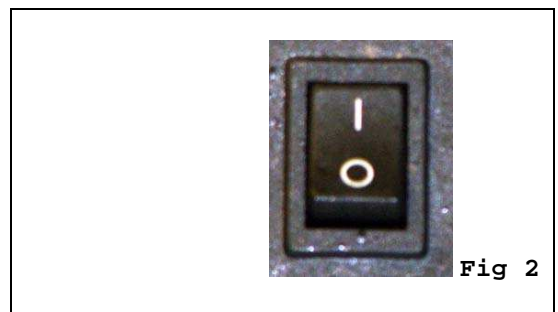


Fig 2



Fig 3